

Technical Data Sheet

EMS Sure Cut – Surfactant

#21630

Epoxy resin blocks are easier to section with glass, as well as, diamond knives when the embedding medium contains 0.5 - 1% of our SURE-CUT. The life of the diamond knife will be extended with its use because it will reduce the friction of the plastic block during sectioning.

Instruction:

Add 0.5- 1% **EMS SURE-CUT** to the embedding resin mixture at the same time of the addition of the curing agent. Mix well at least 20 minutes and embed the specimen as usual.

EMS SURE-CUT works well with all epoxy resin, such as: EMBed-812, Spur Low Viscosity, Araldites etc.