

# Technical Data Sheet

## Technovit® 9100 Histology Technology Accessories

#14655

### Embedding mold specifically for Technovit® 9100




Polyethylene embedding forms 25 mm & "cradle" insert for the round specimen holder on the rotation microtome.

Additional sizes: 15 - 30 - 40 - 50 mm (only cylinder)

#### Application:

1. Remove bottom of form, stick the cradle and the form together

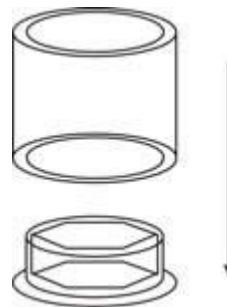

2. Fill up insert




3. Position specimen in the cradle



4. Fill to brim. Turn on vacuum (approx. 10 min in the desiccator) max. 200 mbar, turn off vacuum



5. Put lid on (airtight). Put in the refrigerator at  $-2^{\circ}\text{C}/-4^{\circ}\text{C}$

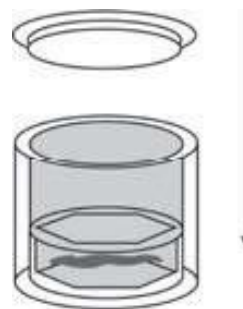

*Source of Information*

Heraeus Kulzer, 2014