# **Johnson's Method For Iron (Microincineration)**

# Catalog #26603

### **Fixation**

10% Buffered Neutral Formalin

# **Sections**

Paraffin at 6 microns

# **Staining**

#### A. Microincineration

- 1. Place unstained paraffin section uncovered on a sheet of asbestos and heat cautiously, using the outer cone from the flame of a Fisher or Meker burner. Keep the flame in motion and heat the slide as uniformly as possible.
- 2. Continue the heating until the specimen, which turns brown then white, has almost completely disappeared. Immediately cover the slide with an inverted pan, (e.g. a pie pan) and allow the slide to cool gradually to rapid cooling may cause the slide to break.
- 3. After the slide has cooled, the completely incinerated specimen is examined under strong oblique lighting with dark field illumination, with or without coverslip.

#### Results

Iron oxide imparts a bright yellow-orange to dark red color to the ash. Specimen must be completely incinerated as carbon particles can also give a yellowish to red color.

### **B. Staining of Incinerated Specimen**

- 1. Dip incinerated specimen in a thin celloidin solution (approx. 0.015% in alcohol: ether, 1:1). Air dry slides for a few minutes and then quickly hydrate to distilled water.
- 2. Stain in working <a href="https://hydrochloric.acid-potassium ferracyanide solution">hydrochloric acid-potassium ferracyanide solution</a>\* for 15 minutes. Rinse in several changes of distilled water. \*Prepare working stain by mixing equal parts of <a href="https://swindle.com/5%/bydrochloric.acid">5% bydrochloric acid</a> and <a href="https://swindle.com/5%/bydrochloric.acid">5% potassium ferracyanide</a> just before use.
- 3. Air dry or dry in a 60°C oven.
- 4. Mount with Permount.

## C. Direct observation of Prussian Blue Reaction on the Ash

- 1. Create a well paraffin around the specimen.
- 2. Pour working <u>Hydrochloric Acid-Potassium Ferracyanide Solution</u> (for preparation, sheet step B-2) into the paraffin well and observe under the microscope.

# Staining Results (B & C)

Iron Oxide	Blue
------------	------

# References

Fenton, R.H., Johnson, F.B., and Zimmerman, L.E., J. Histo-chem. Cytochem., 12:153:155, 1964.

L una, L.G., (ed.), Manual of Histologic Staining Methods of the AFIP , 3 rd edition, McGraw-Hill, NY., p. 181., c 1968.